



Mounting Procedure of FO Air Exhausting Device


1. Take out the ceramic plug of exhaust port at the upside of equipment.


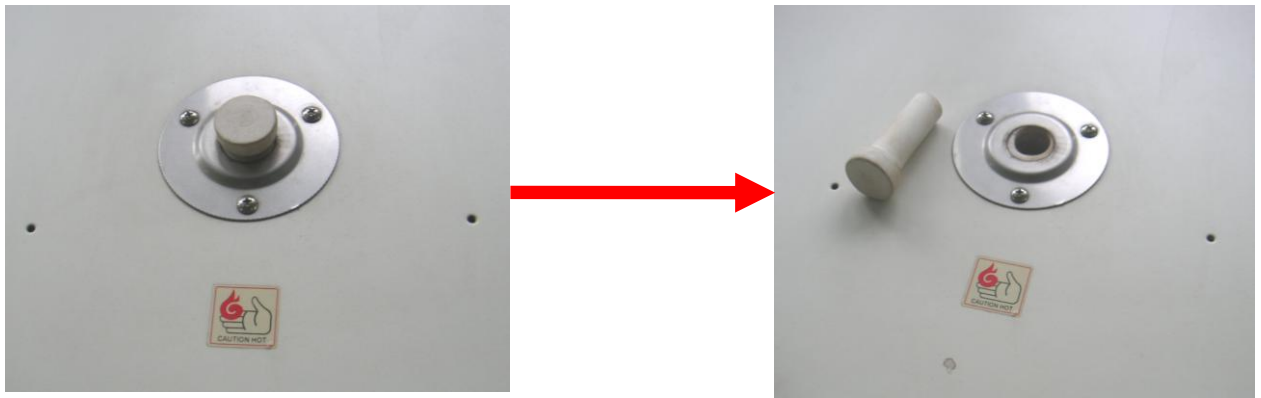


2. Put the air exhausting device on the equipment, corresponding to threaded holes.

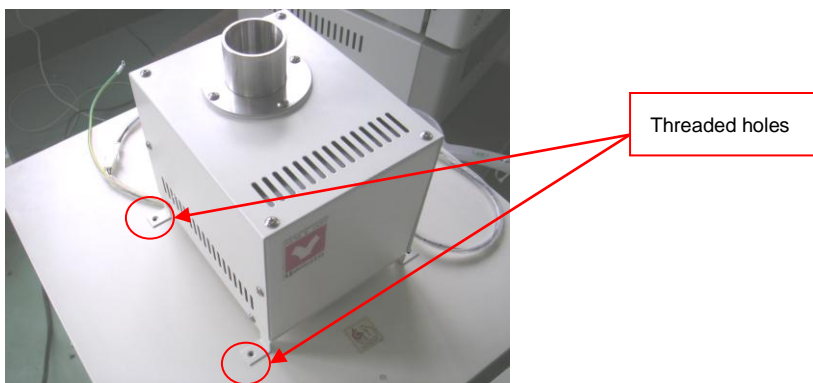


3. Fix the air exhausting device.

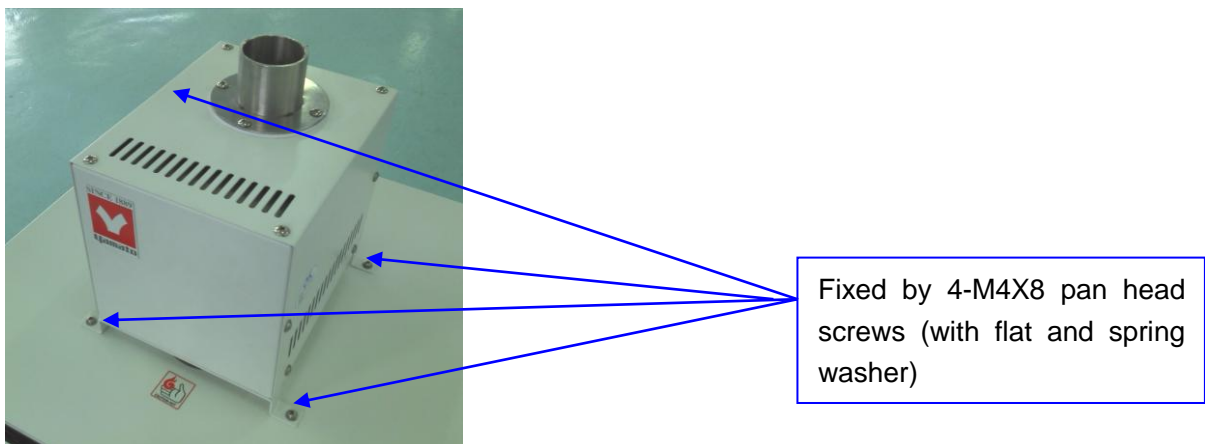


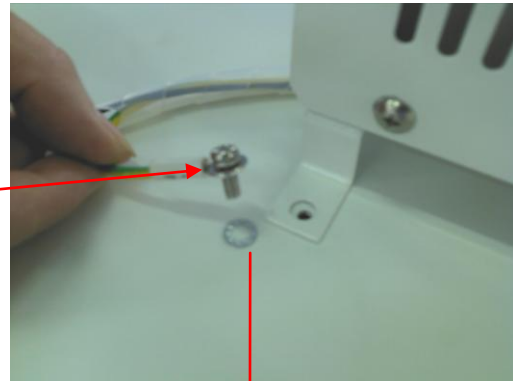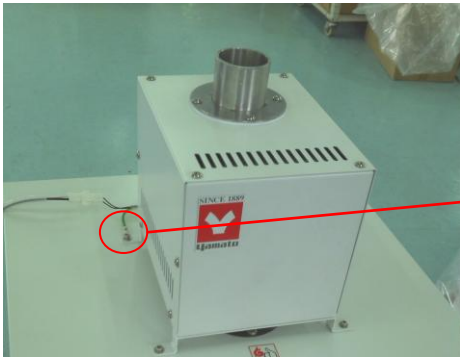


Before fixing left-rear screw, put internal tooth washer to ground between the main unit and air exhausting device, and then tighten the screw.

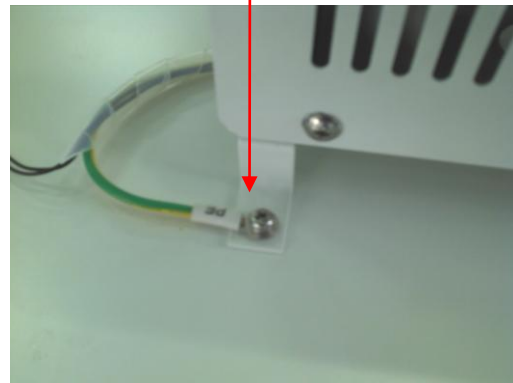

4. Wiring

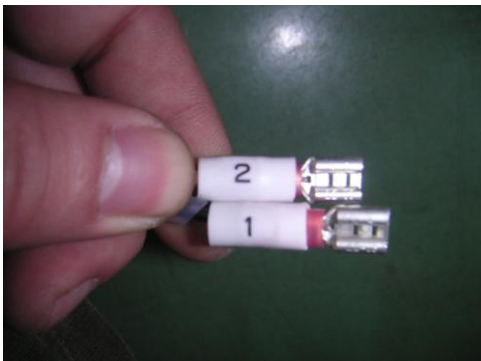

- Disassemble the underside electric backboard;
- According to the wire marking of accessory wire harnesses, insert into corresponding terminal blocks;
- Confirm the wiring is correct and firm, and then mount the electric backboard.

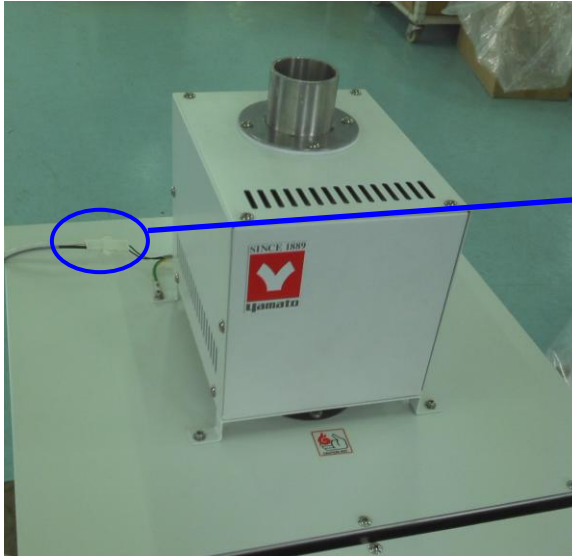

Connect the outgoing lines of equipment and air exhausting device.



Disassemble the original screws, and use accessory 2-M4X8 big flat head screws to fix the cable clips. Do not screw tightly too much.



5. Complete assembly.



Lead out the wire harness of
air exhausting device here

Notice: when completing assembly, if the equipment is power-on, the air exhausting device will run at once. If not use the air exhausting device, please pull out the wiring of step 4, and take the exhausting device away from the equipment.

Parts detail list

NO.	Name	Qty.
1	Air exhausting device	1
2	Cross-head double washers screw M4X8	4
3	Internal tooth washer $\Phi 4$	1
4	Big flat head screw M4X8 (attached when separately delivering the air exhausting device)	2
5	Cable clip (attached when separately delivering the air exhausting device)	2
6	Wire harness (attached when separately delivering the air exhausting device)	1
7	Mounting procedure	1

Notice: Part 4, 5 and 6 have already been assembled at FO main unit, and the big flat head screw, cable clip and wire harness are attached only when separately delivering the air exhausting device.