


SN / SQ QUICK OPERATION GUIDE

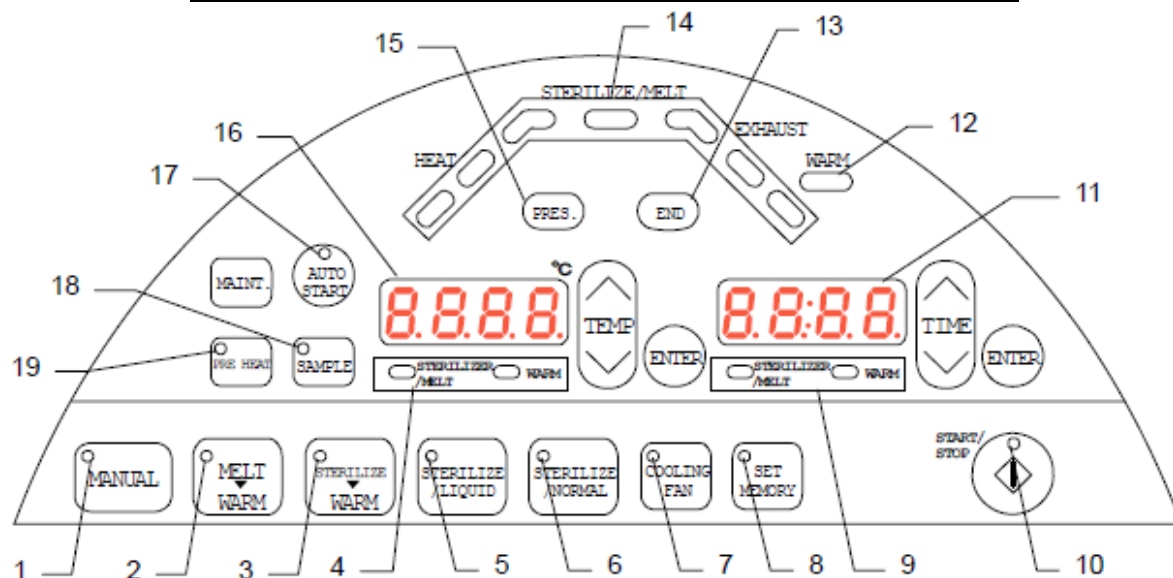

No.	Name	Function
1	Manual course lamp (MANUAL)	This lamp lights up when manual course is selected.
2	Melting & Warm course lamp (MELT→WARM)	This lamp lights up when melting & warm course is selected.
3	Sterilization & Warm course lamp (STERILIZE→WARM)	This lamp lights up when sterilization & warm course is selected.
4	Sterilization/melting & warm (STERILIZER/MELT, WARM) temperature lamp	This lamp blinks at temperature setting of sterilization/melting & warm.
5	Liquid sterilization course lamp (STERILIZE/LIQUID)	This lamp lights up when liquid sterilization course is selected.
6	Apparatus sterilization course lamp (STERILIZE/NORMAL)	This lamp lights up when apparatus sterilization course is selected.
7	Forced cooling lamp (COOLING FAN)	This lamp lights up when forced cooling process is selected.
8	MEMORY lamp (SET MEMORY)	This lamp lights up at the setting of memory.
9	Sterilization/melting & warm (STERILIZER/MELT, WARM) time lamp	This lamp blinks at the setting of time for sterilization/melting & warm
10	Start/stop lamp (START/STOP)	This lamp lights up during operation and goes out at operation stop.
11	Time display screen	The display blinks at the setting of time parameters. It displays the remaining time during operation of equipment.
12	Heat-retention lamp (WARM)	This lamp blinks at the setting of heat-retention process.
13	End lamp (END)	This lamp blinks when operation is completed.
14	Operation monitoring lamp	Refer to Page 20.
15	Pressurization lamp (PRES.)	This lamp lights up when the equipment is being pressurized.
16	Temperature display screen	The display blinks at the setting of temperature parameters. It displays the temperature inside the chamber during operation of equipment.
17	Auto start lamp (AUTO START)	This lamp lights up when auto start key is pressed. It blinks after the auto start time is set.
18	Sample temperature lamp (SAMPLE)	This lamp lights up when the sample temperature key (SAMPLE) is selected when optional sample temperature sensor, which is sold separately, is attached.
19	Preheating lamp (PRE HEAT)	This lamp lights up when the preheating key (PRE HEAT) is selected. It blinks during the preheating operation.