


## HIGH ALTITUDE STERILIZER ADJUSTMENT

### I. Modify Steam Saturation Temp

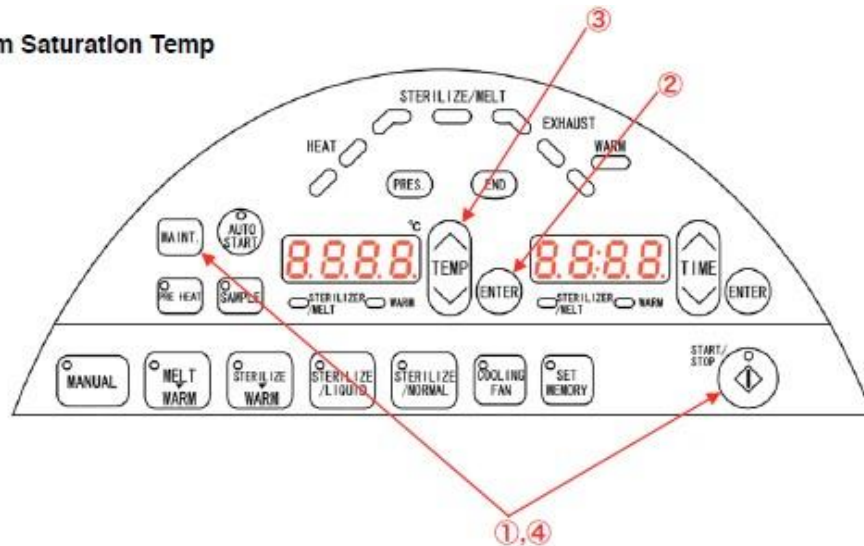

- ① Press "MAINT", then "START/STOP" and hold them down together for 2 seconds (enter factory setting mode).
- ② Press "ENTER" repeatedly until "SAtE" can be viewed in the TIME display.\*
- ③ Enter a temperature value of 89 into the TEMP display, using "TEMP" arrow keys. If value is already 89, leave it as is.

\*If setting "SAtE" cannot be found, follow the "unhide settings procedure" below.

### II. Modify Steam Saturation Temp Range

- ① Skip to next step (stay in factory setting mode)
- ② Press "ENTER" repeatedly until "SAtb" can be viewed in the TIME display.\*
- ③ Enter a range value of 3.0, using "TEMP" arrow keys. If value is already 3.0, leave it as is.

\*If setting "SAtb" cannot be found, follow the "unhide settings procedure" below.

### III. Modify Steam Saturation Temp Time

- ① Skip to next step (stay in factory setting mode)
- ② Press "ENTER" repeatedly until "SAt i" can be viewed in the TIME display.\*
- ③ Enter a time value of 1, using "TEMP" arrow keys. If value is already 1, leave it as is.
- ④ Press "MAINT" to return to normal operation mode.

\*If setting "SAt i" cannot be found, follow the "unhide settings procedure" below.

### Unhide Settings Procedure

- ① Press "MAINT", then "START/STOP" and hold them down together for 2 seconds **while in factory setting mode** (during steps I, II, or III above).
- ② Press "ENTER" repeatedly until the desired setting ("SAtE", "SAtb" or "SAt i") can be viewed in the time (right) display.
- ③ Choose "oFF" using the "TEMP" arrow key.
- ④ Press and hold "MAINT" and "START/STOP" together for 2 seconds again to return to factory setting mode.