



**Glass Set
for
Rotary Evaporator**

**RG202A-A
RG202B-A**

Instruction Manual

First Edition

Yamato Scientific Co., Ltd.

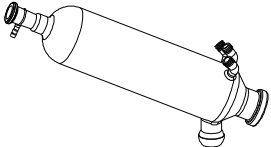
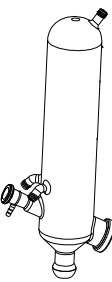
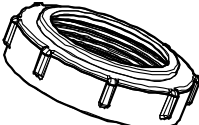


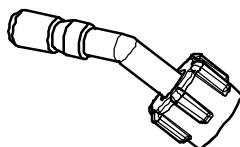
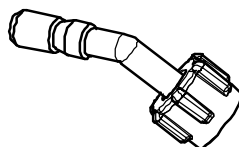
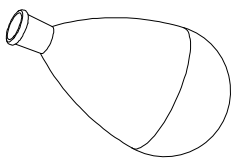
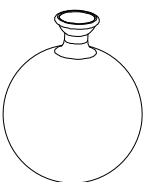
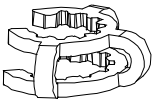
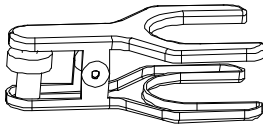
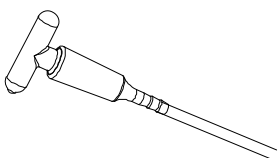
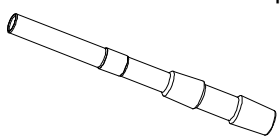
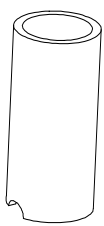

Printed on recycled paper

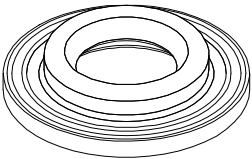
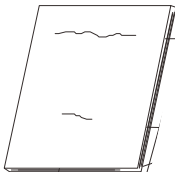
1. Compatible Models

Rotary Evaporator RE202-A/212-A

2. Packing List

Check before operation that all the accessories are complete. Contact original dealer of purchase if anything is missing.

<p>①-A Cooling condenser type A 1 pc</p>  <p>* RG202A-A only</p>	<p>①-B Cooling condenser type B 1 pc</p>  <p>* RG202B-A only</p>		<p>②Condenser nut 1 pc</p> 
<p>③Coil ring 1 pc</p> 	<p>④Flask removal tool 1 pc</p>  <p>Standard: ¥24</p>	<p>⑤Vacuum nozzle (gray) 1 set</p> 	<p>⑥Cooling nozzle (black) 2 set</p> 
<p>⑦Evaporation flask 1000 mL 1pc</p>  <p>Standard: ¥24/40</p>	<p>⑧Receiving flask 1000 mL 1pc</p>  <p>Standard: S35/20</p>	<p>⑨Flask clip 1 pc</p>  <p>¥24</p>	<p>⑩Ball joint clamp 1 pc</p> 
<p>⑪Sample feed stopcock 1 pc</p>  <p>Standard: ¥19/38 * PTFE tube included</p>	<p>⑫Rotary joint 1 pc</p>  <p>RG202A-A : L286 mm RG202B-A : L210 mm Standard: ¥24/40</p>	<p>⑬Condenser insulation 1 pc</p> 	<p>⑭Nozzle insulation 2 pcs</p> 

<p>⑮ Vacuum seal</p> <p>1 pc</p>  <p>Material: FKM</p>	<p>⑯ Instruction manual</p> <p>1 copy</p> 
---	---

* Carefully read RE202-A/212-A standard instruction manual, and set up glass set properly for safe operation.



Use caution when handling glassware

Any damage on glassware may lead to serious accidents.
Check glassware for damage before operation.

Limited Liability

Always operate equipment in strict compliance to the handling and operation procedures set forth by this instruction manual.

Yamato Scientific Co., Ltd. assumes no responsibility for malfunction, damage, injury or death, resulting from negligent equipment use.

Never attempt to disassemble, repair or perform any procedure on RE units which are not expressly mandated by this manual. Doing so may result in equipment malfunction, serious personal injury or death.

Notice

- Instruction manual descriptions and specifications are subject to change without notice.
- Yamato Scientific Co., Ltd. will replace flawed instruction manuals (pages missing, pages out of order, etc.) upon request.

Instruction Manual

Glass Set for Rotary Evaporator

RG202A-A

RG202B-A

First Edition: July 5, 2019

Revised:

Manufacturer

Yamato Scientific Co., Ltd.

Harumi Triton Square Office Tower Y (36F)

1-8-11 Harumi, Chuo-ku, Tokyo 104-6136, JAPAN

Yamato Scientific America Inc.

925 Walsh Avenue, Santa Clara,

CA 95050, U.S.A

<http://www.yamato-usa.com>

Toll Free: 1-800-2-YAMATO (1-800-292-6286)